


Simple Concentration and Analysis of Fragrance Components in Dairy Products

- Using Sorptive Media MonoTrap Sampler

MonoTrap RCC18 (with activated carbon) was used to collect flavor components from dairy products using a head-space method (HS).


Preliminary processing procedure

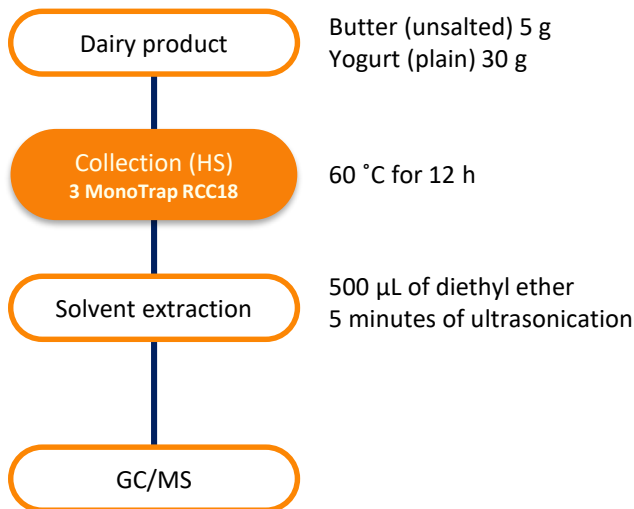

GC Conditions

System : GC - MS
Column : InertCap Pure-WAX
 0.25 mm I.D. x 30 m df = 0.25 µm
Col. Temp. : 40 °C (5 min) - 4 °C/min - 220 °C
Carrier Gas : He 1 mL/min
Injection : Splitless 250 °C
Detection : MS Scan (35-500 m/z)
Sample Size : 1 µL


Yogurt

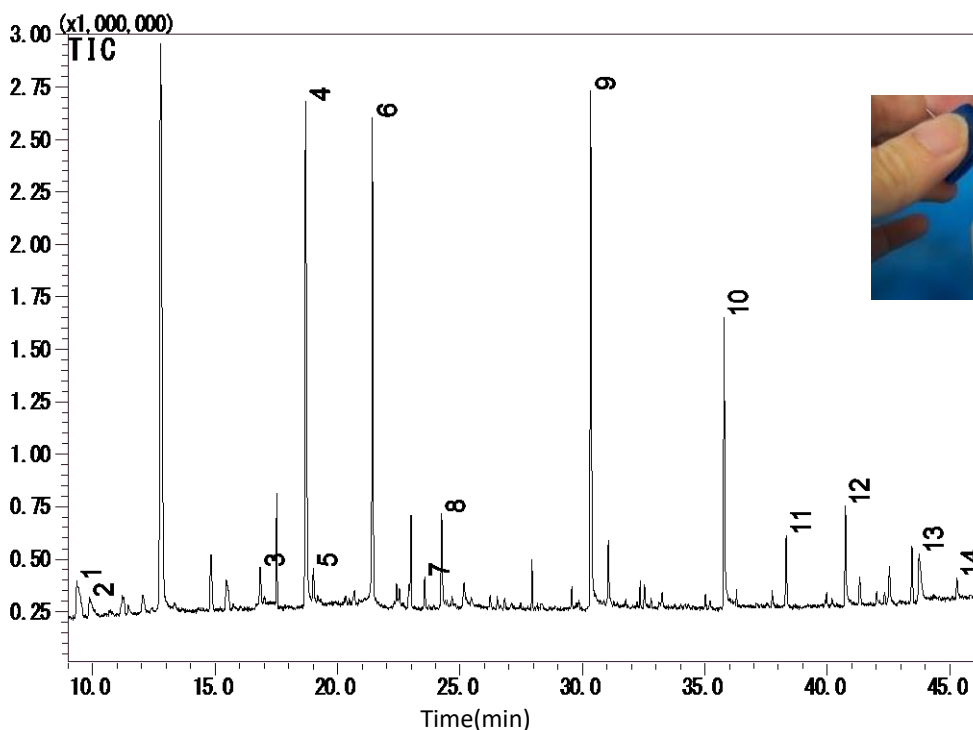

- | | |
|------------------------------------|----------------------------|
| 1. 2-Heptanone | 8. Butanoic acid |
| 2. D-Limonene | 9. Hexanoic acid |
| 3. 2-Nonanone | 10. Octanoic acid |
| 4. Acetic acid | 11. δ-Decalactone |
| 5. Furfural | 12. N-Decanoic acid |
| 6. (3-Methyl-oxiran-2-yl)-methanol | 13. Benzenecarboxylic acid |
| 7. 2-Undecanone | 14. Dodecanoic acid |

*According to library search results
Red letters; identification results with standard samples

Butter

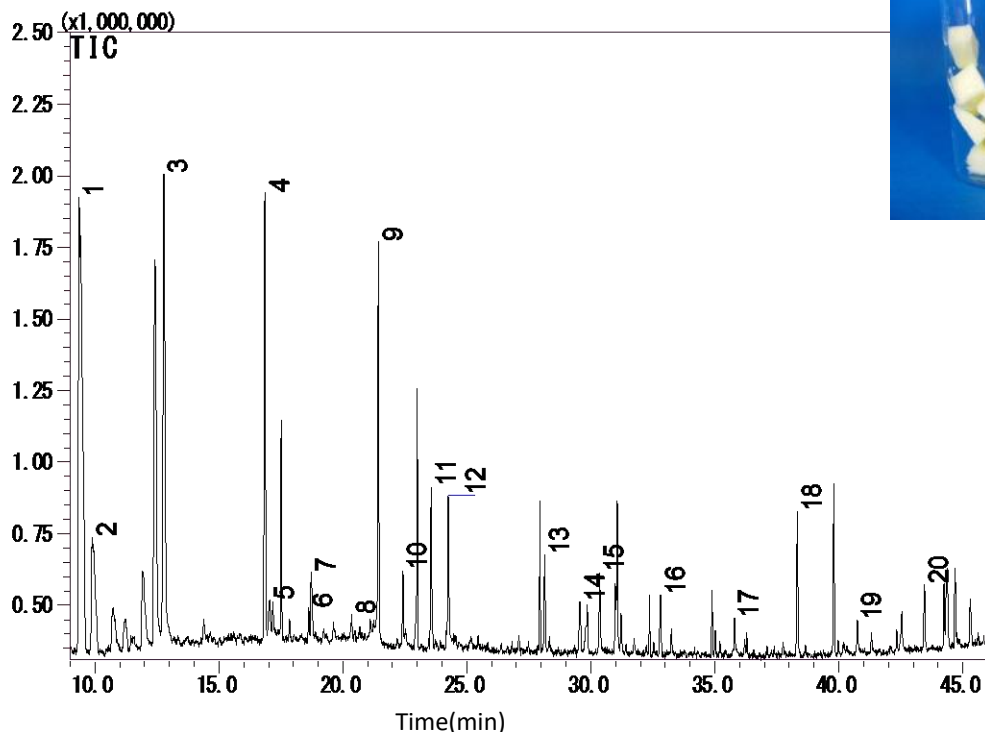

- | | |
|---------------------------|---------------------------------|
| 1. 2-Heptanone | 11. 2-Undecanone |
| 2. D-Limonene | 12. Butanoic acid |
| 3. Unknown | 13. δ -Caprolactone |
| 4. 2-Nonanone | 14. 2-Tridecanone |
| 5. Nonanal | 15. Hexanoic acid |
| 6. Ethyl caprylate | 16. δ -Octalactone |
| 7. Acetic acid | 17. Octanoic acid |
| 8. Ethylhexanol | 18. δ -Amylvalerolactone |
| 9. 1,3-Butanediol | 19. N-Decanoic acid |
| 10. 1-Octanol 2-Heptanone | 20. δ -Dodecalactone |

*According to library search results

Red letters; identification results with standard samples

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome
Shinjuku-ku, Tokyo,
163-1130, Japan
Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences B.V.

De Sleutel 9
5652 AS Eindhoven
The Netherlands
Phone: +31 (0)40 254 95 31
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
Phone: 310-265-4424
Fax: 310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,
Far East International Plaza,
NO,317 Xianxia Road,
Changning District.
Shanghai, China P.C. 200032
Phone: +86 (0)21-6278-2272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn

International Distributors

Visit our Website at www.glsciences.com/distributors