


Analysis of Paroxetine hydrochloride

(Under the Condition of the Japanese Pharmacopoeia)

Data No. LB206-0894

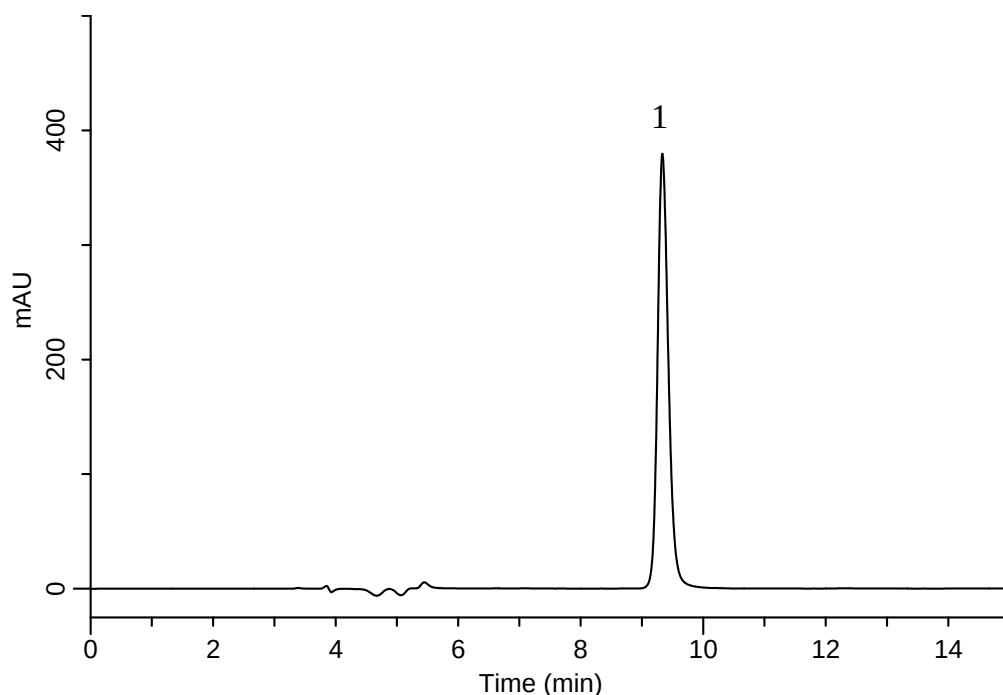

Conditions

System : GL-7400 HPLC system
Column : Inertsil TMS
(5 μ m, 250 x 4.6 mm I.D.)
Column Cat. No. : 5020-18161
Eluent : A) CH₃CN
B) Buffer*
C) Triethylamine
A/B/C = 40/60/1, v/v/v (pH5.5, CH₃COOH)
Flow Rate : 0.7 mL/min
Col. Temp. : 30 °C
Detection : UV 295 nm (GL-7452 PDA Detector)
Injection Vol. : 10 μ L
Sample : Standard

*Dissolve 3.85 g of ammonium acetate in 1000 mL of water.
Adjust pH 4.5 by acetic acid.

Analyte:

1. Paroxetine Hydrochloride 500 mg/L

Theoretical plate number : 14,356 ($\geq$ 5,000)
Tailing factor : 1.20 ($\leq$ 2.0)
RSD of the peak area (%) (n=6) : 0.36 ($\leq$ 1.0)