


Residual solvents in pharmaceuticals Class1

Data No. GA089-0000

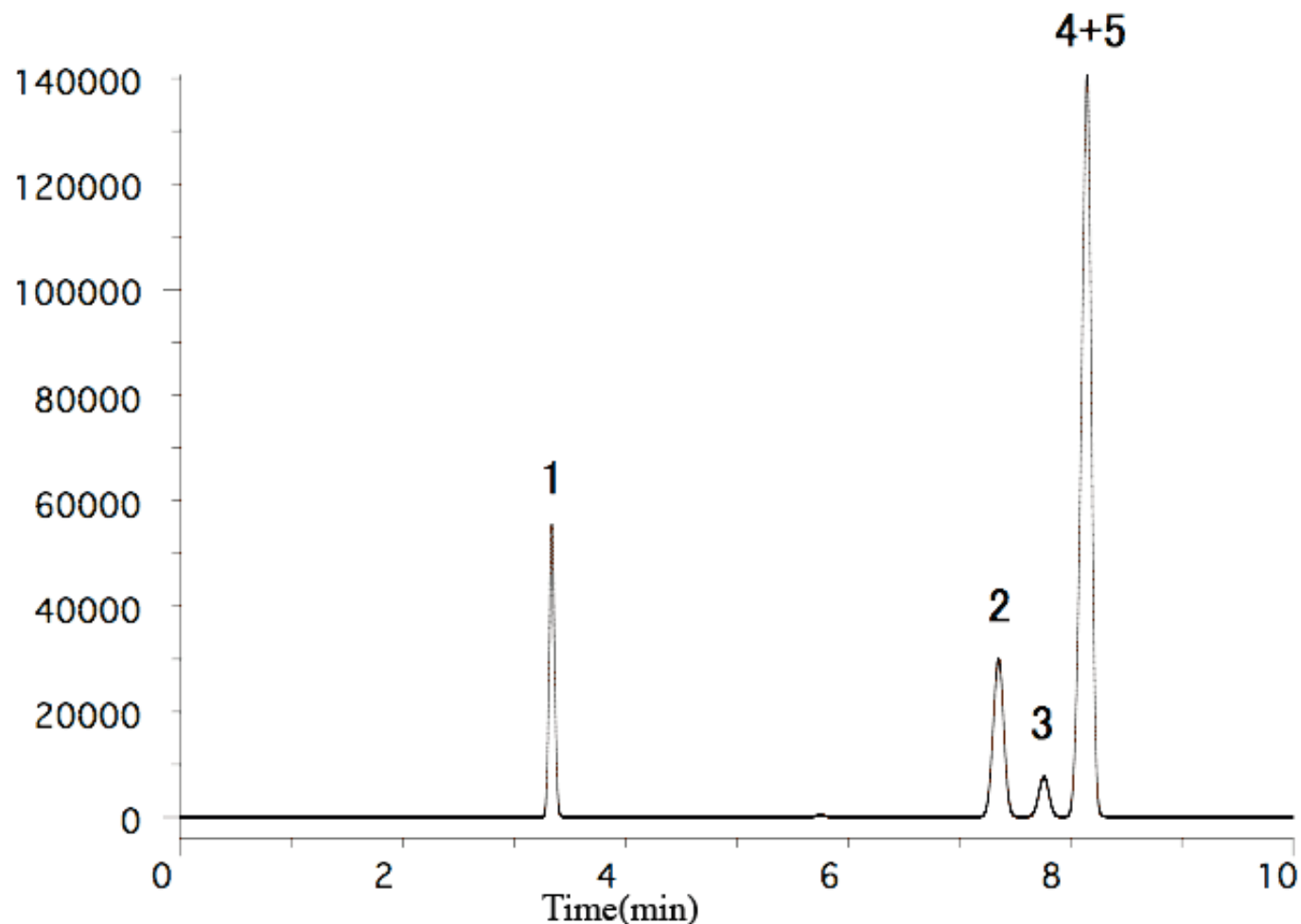

System : GC/FID
Column : InertCap624
0.53mm I.D. × 30m df=3µm
Column Temp. : 40°C (5min) – 5°C/min – 65°C
Carrier Gas : He 20kPa 35cm/sec
Injection : Split Flow 200ml/min
240°C
Detection : FID Range 10²
240°C
Sample : Solvents
Mixed evenly
Sample Size : 0.4µL

Solutes 1. 1,1-Dichloroethylene
2. 1,1,1-Trichloroethane
3. Carbon tetrachloride
4. 1,2-Dichloroethane
5. Benzene