EXHAUST VALUES • LUFTMÄNGDER

Sugearme / [Sugearme](#)

Ø 50 mm	Normal	65 m ³ /h	Minimum	45 m ³ /h	Maksimal	85 m ³ /h
Ø 63 mm	Normal	90 m ³ /h	Minimum	60 m ³ /h	Maksimal	120 m ³ /h
Ø 75 mm	Normal	140 m ³ /h	Minimum	80 m ³ /h	Maksimal	180 m ³ /h
Ø100 mm	Normal	270 m ³ /h	Minimum	140 m ³ /h	Maksimal	400 m ³ /h

Fixed airflow / [Fast luftstrøm](#)

Sugeskabe / [Sugeskabe](#)

LAV / BAB	Width / Bredde	900 mm	Normal	400 m ³ /h
LAV / BAB	Width / Bredde	1200 mm	Normal	500 m ³ /h
LAV / BAB	Width / Bredde	1500 mm	Normal	600 m ³ /h
LAV / BAB	Width / Bredde	1800 mm	Normal	800 m ³ /h

Fixed airflow / [Fast luftstrøm](#)

With VAV system, see fume cupboards / [Med VAV system tilsluttet, se under stinkske](#)

Chemical & Toxic cabinets / **Kemikalie & Giftskabe**

There is typically a requirement that the air in a chemical cabinet must be replaced at a certain frequency to prevent the accumulation of hazardous fumes. A common recommendation is an air course of at least 10 times per hour.

There must be continuous ventilation to ensure constant removal of potentially hazardous fumes from the chemical cabinet. This can either be an independent ventilation system for each cabinet or a central system that serves multiple cabinets.

If flammable or explosive chemicals are stored, the ventilation must be explosion-proof. When storing corrosive substances, the ventilation system must be corrosion resistant.

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Normal 160 m3/h
Normal 200 m3/h
Normal 300 m3/h

Fireproof Chemical cabinets / Brandsikre kemikalieskabe

For fireproof cabinets, especially those that meet the EN 14470-1 standard, it is recommended that cabinets with flammable liquids have an active ventilation system. This system must ensure a constant exchange of air inside the cabinet, which is typically at least 10 times per hour. This corresponds to a ventilation rate of approx. 1-2 m³/h, depending on the size and contents of the cabinet.

If the cabinet stores flammable or explosive chemicals, it is important that the ventilation system is explosion-proof to prevent ignition of the fumes.

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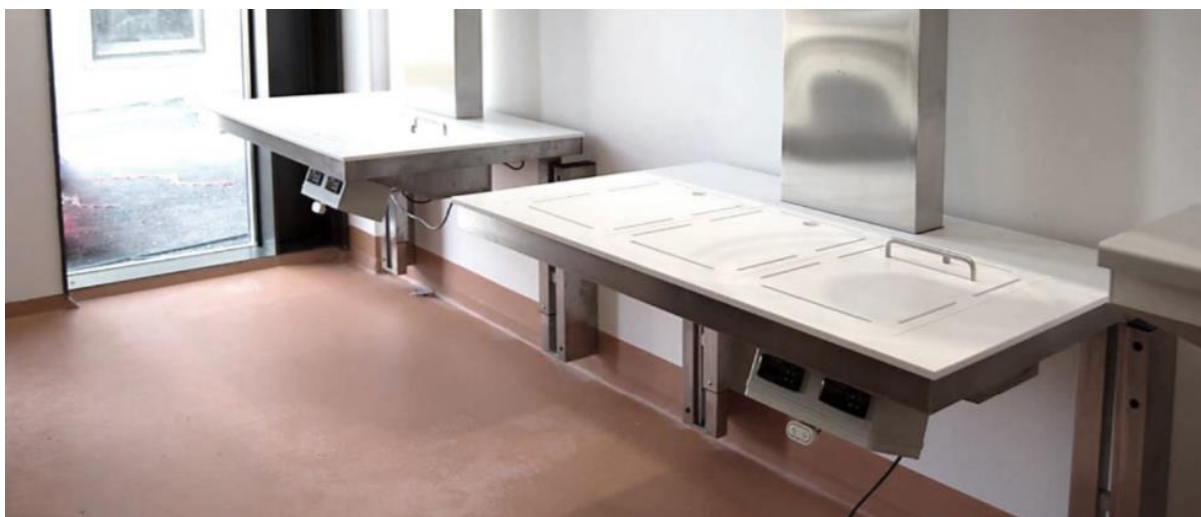
Exhaust: **300 m³/h**

Pressure drop / [Tryktab](#): 9 Pascal

Safe outback from work area up to 4 cm over the ventilated area

Exhaust: **500 m³/h**

Pressure drop / [Tryktab](#): 15 Pascal

Safe outback from work area up to 8 cm over the ventilated area

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LABMODUL A/S