

# Concentration analysis of toothpaste flavor components - Using MonoTrap Simple Enrichment Tools

MonoTrap RGC18TD, a simple enrichment tool was used to screen and analyze the flavor components in toothpaste. MonoTrap RGC18TD is dedicated to heat desorption sampling, it is supplied individually packaged in ampoules and does not require conditioning prior to use.

## Preliminary processing procedure

Toothpaste

Put 1 g (approximately one dose) into a 40 mL vial.


Collection (HS)  
One MonoTrap RGC18TD

36 °C for 10 minutes

TD/GC/MS

**Conditions**

<b>System</b>	: GC – MS - Thermal Desorption
<b>Column</b>	: InertCap Pure-WAX ProGuard 0.32 mm I.D. x 2 m + 120 m, df = 0.50 µm
<b>Col.Temp.</b>	: 40 °C (3.5 min) - 4 °C/min - 250 °C (20 min)
<b>Carrier Gas</b>	: He 2.5 mL/min (constant flow)
<b>Desorb Temp.</b>	: 200 °C
<b>Time</b>	: 5 min
<b>Mode</b>	: Split (Split ratio 5:1)
<b>Cryo Trapping</b>	: -150 °C
<b>Injection Temp.</b>	: 250 °C
<b>Detection</b>	: MS Scan ( <i>m/z</i> : 29 - 800)

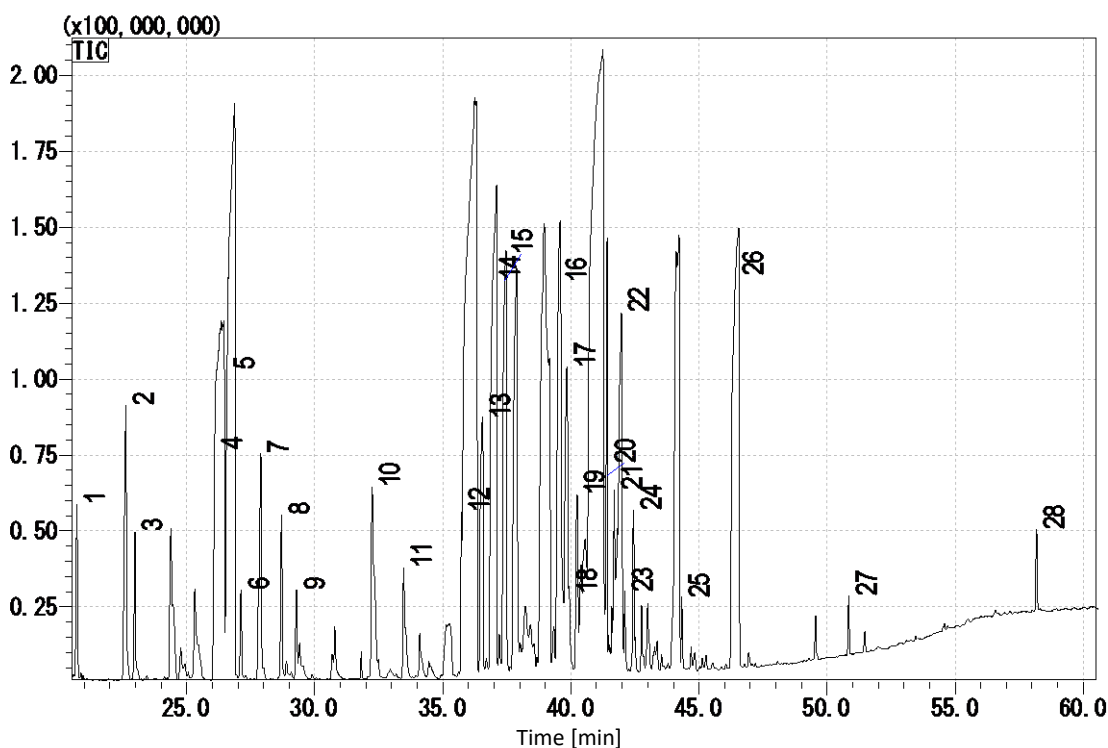

1	Camphene	11	Camphor	21	Isomenthol
2	Fenchene	12	Ethyl hexanol	22	Estragole
3	Sabinen	13	Cineron	23	Dihydrocarvyl acetate
4	Limonen	14	Menthone	24	Terpineol
5	Cineol	15	Dihydrocarveol	25	Anethol
6	Ocimene	16	Menthol	26	Estragole
7	Terpinolen	17	Terpineol	27	Menthyl lactate
8	Menthatriene	18	Bornyl acetate	28	Thymol
9	Terpinolene	19	Isopulegol		
10	Octanol	20	Caryophyllene		

\* According to library search results

## Products

### Absorbent

Description : MonoTrap RGC18TD

Cat.No. : 1050-74201

\* Shipped in ample tube.



MonoTrap RGC18TD



MonoTrap Trial Kit for TD

Contents	Cat.No.
<ul style="list-style-type: none"> <li>● MonoTrap TD 10 pcs (Select one from 3 types as below) MonoTrap RSC18 TD, RGC18 TD, RGPS TD</li> <li>● Glass Tube 1 pc ( Select a tube for OPTIC / LINE, or for T-Dex / ATD / TD-20)</li> <li>● MT Holder 1 pc</li> <li>● CLEAN PIN HOLE SEPTUM WITH VIAL 1 pc</li> </ul>	1050-07800

### GC Capillary Columns

Description : InertCap Pure-WAX

Size : 0.32 mm I.D. x 2 m (ProGuard) + 120 m, df = 0.50  $\mu$ m

Contact us for detail



InertCap Pure-WAX

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

#### GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome  
Shinjuku-ku, Tokyo,  
163-1130, Japan  
Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsciences.co.jp](mailto:world@glsciences.co.jp)  
Web: [www.glsciences.com](http://www.glsciences.com)

#### GL Sciences B.V.

De Sleutel 9  
5652 AS Eindhoven  
The Netherlands  
Phone: +31 (0)40 254 95 31  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

#### GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
Phone: 310-265-4424  
Fax: 310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

#### GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,  
Far East International Plaza,  
NO,317 Xianxia Road,  
Changning District.  
Shanghai, China P.C. 200032  
Phone: +86 (0)21-6278-2272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)

#### International Distributors

Visit our Website at [www.glsciences.com/distributors](http://www.glsciences.com/distributors)