


# Easy Enrichment of VOC from Papers Before & After Printing with MonoTrap

MonoTrap is a hybrid novel adsorbent that combines large surface area and properties of silica gel, activated carbon, and ODS. Utilizing the surface of porous silica, it can be used as a tool that makes use of the large surface area and the adsorbent effect caused by the inclusion of activated carbon for high collection efficiency and for short time and sensitive analysis. MonoTrap DCC18 (with activated carbon) was used to analyze VOCs in new vehicles. Activated carbon is effective in collecting polar and low boiling components. It is also easy to collect by simply placing a small collection material inside the car room. Capture concentration can also be used as usual, and simple and stress-free enrichment analysis can be performed.




## Protocol

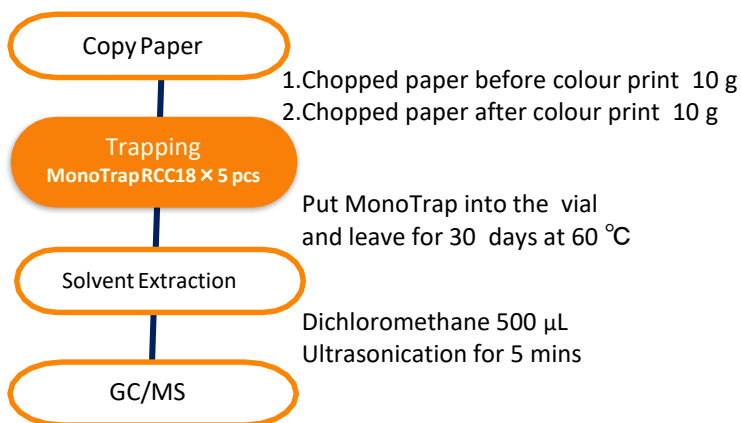

### GC Conditions

<b>System</b>	: GC/MS
<b>Column</b>	: InertCap Pure-WAX 0.25 mm I.D. × 30 m df = 0.25 µm
<b>Col. Temp.</b>	: 40 °C (5 min) - 4 °C /min - 250 °C
<b>Carrier Gas</b>	: He 1 mL/min
<b>Injection</b>	: Splitless 250 °C
<b>Detection</b>	: MS Scan ( <i>m/z</i> 35-500)
<b>Sample Size</b>	: 1 µL

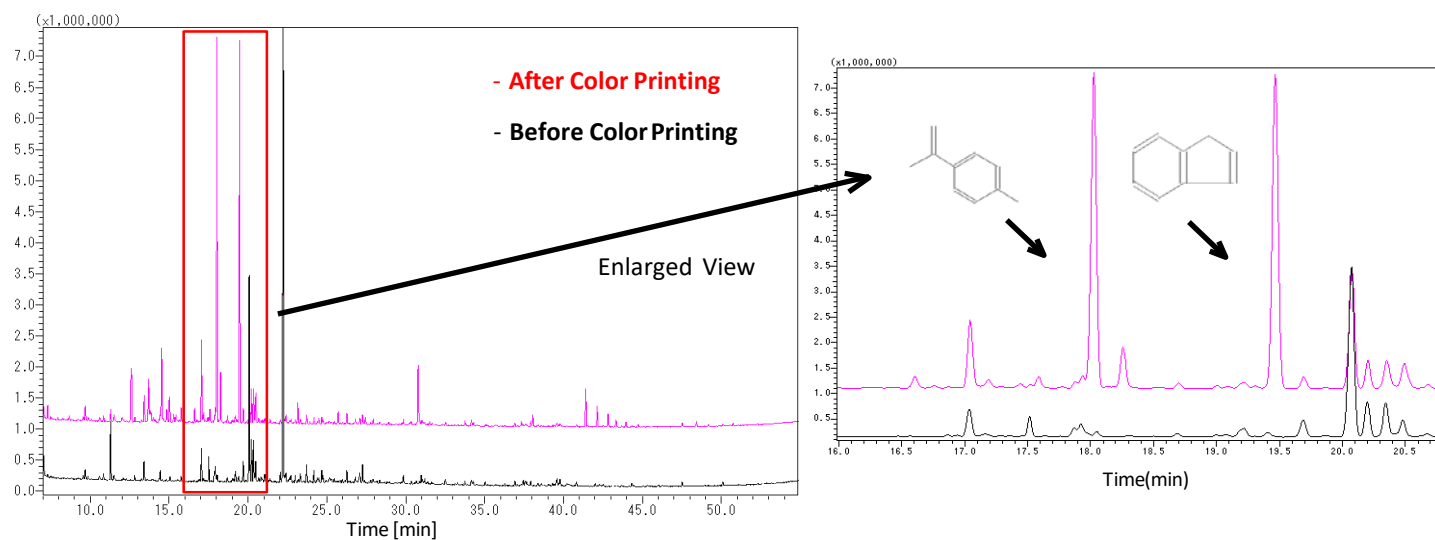

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

### GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome  
Shinjuku-ku, Tokyo,  
163-1130, Japan  
Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsc.co.jp](mailto:world@glsc.co.jp)  
Web: [www.glsciences.com](http://www.glsciences.com)

### GL Sciences B.V.

De Sleutel 9  
5652 AS Eindhoven  
The Netherlands  
Phone: +31 (0)40 254 95 31  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

### GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,  
Far East International Plaza,  
NO,317 Xianxia Road,  
Changning District.  
Shanghai, China P.C. 200051  
Phone: +86 (0)21-6278-2272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)

### GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
Phone: 310-265-4424  
Fax: 310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

### International Distributors

Visit our Website at:

<https://www.glsciences.com/company/distributor.html>