


Simple Concentration and Analysis of Spice Flavor Components - Using Sorptive Media MonoTrap

In this application note, MonoTrap RCC18 (with activated carbon) was used to simply concentrate and analyze the flavor components of spices.

Preliminary processing procedure

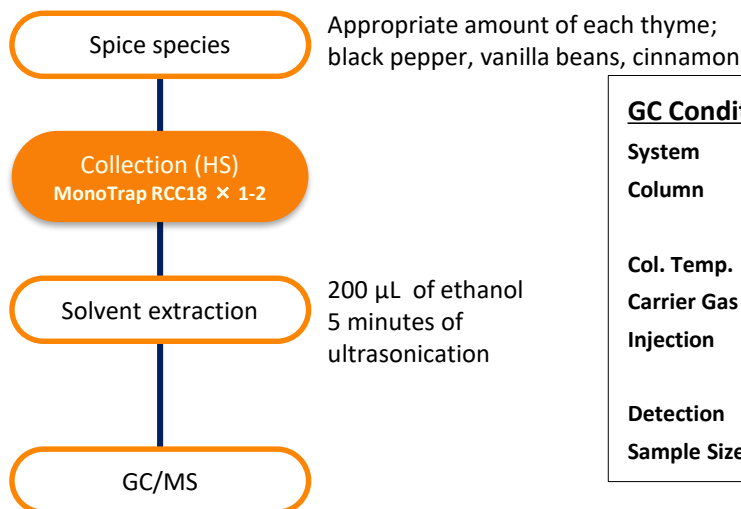


GC Conditions

System : GC - MS
Column : InertCap Pure-WAX
 0.25 mm I.D. x 30 m df = 0.25 µm
Col. Temp. : 40 °C(5 min) - 5 °C/min - 250 °C
Carrier Gas : He 1 mL/min
Injection : Split 1:20
 250 °C
Detection : MS Scan (m/z : 35-600)
Sample Size: 1 µL



Image of collection vials

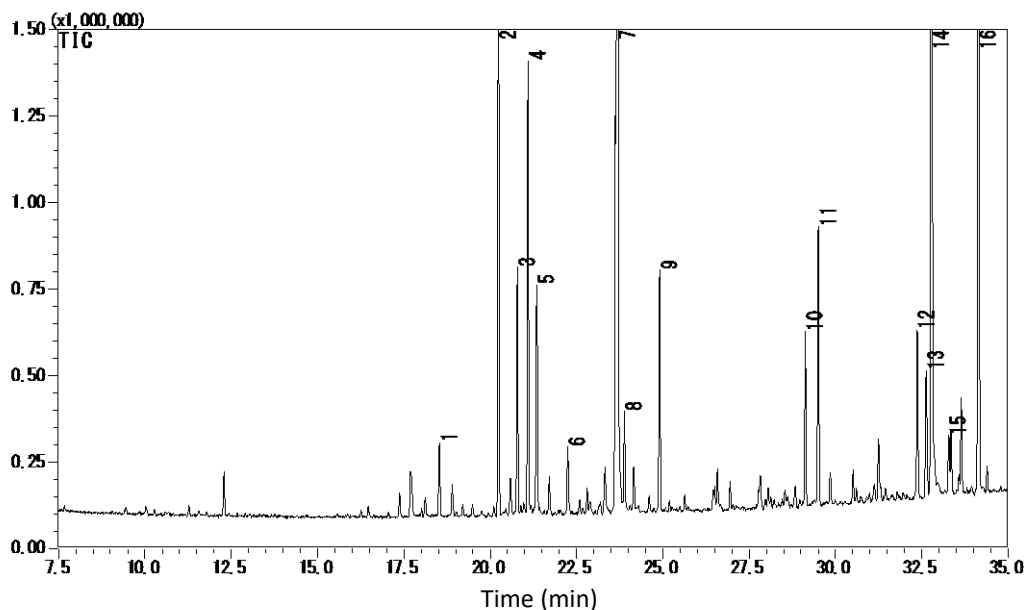


1. Copaene
2. β -Linalool
3. 1,7,7-Trimethylbicyclo[2.2.1]hept-2-yl acetate
4. 4,11,11-Trimethyl-8-methylenebicyclo[7.2.0]undec-4-ene
5. *p*-Menth-1-en-4-ol
6. Alloaromadendrene
7. Borneol
8. 4,7,7-Trimethylbicyclo[4.1.0]heptan-3-ol

9. γ -Muurolene
10. 3,7-Octadiene-2,6-diol, 2,6-dimethyl-
11. Caryophyllene oxide
12. Spathulenol
13. 2,6-Dimethyl-1,7-octadiene-3,6-diol
14. Unknown
15. Tau.-Cadinol
16. Carvacrol

*According to the library search results


Black pepper

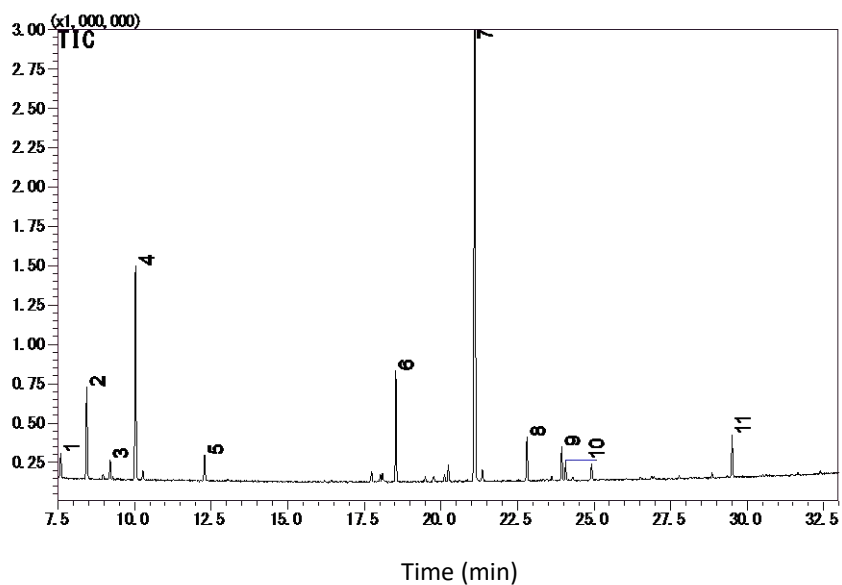

1. 4(10)-Thujene
2. 3-Carene
3. β -Myrcene
4. D-Limonene
5. *o*-Cymene
6. Copaene
7. Caryophyllene
8. α -Caryophyllene
9. Eudesma-4(14),11-diene
10. α -Selinene
11. Caryophyllene oxide

*According to library search results


Vanilla beans

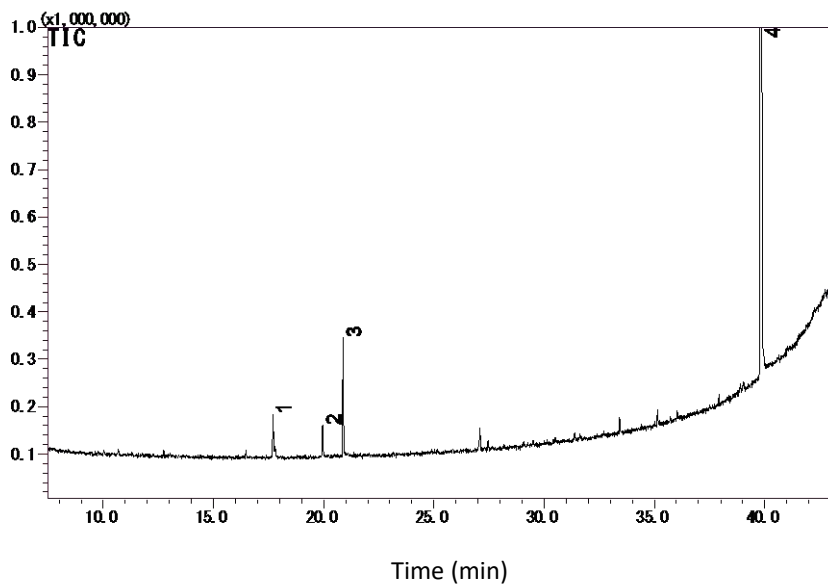

1. Acetic acid
2. Butanediol
3. Butanediol
4. Vanillin

*According to library search results

Cinnamon

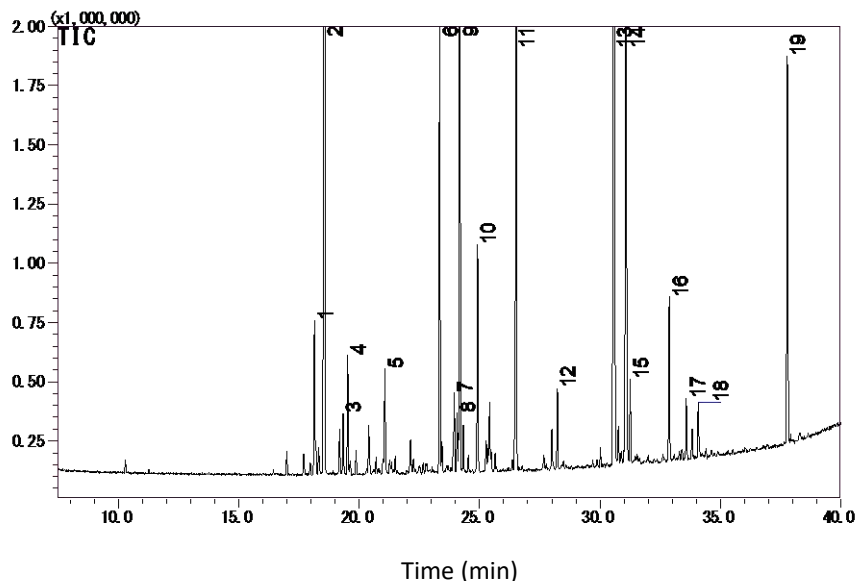

1. Cycloisosativene
2. α -Cubebene
3. Sativen
4. Sativen
5. β -Elemene
6. γ -Muurolene
7. Eudesma-4(14),11-diene
8. β -Chamigrene
9. α -Muurolene
10. δ -Cadinene
11. Calamenene
12. α -Calacorene
13. Cinnamaldehyde
14. 3-Methyl-7,8-dihydroquinolin-5(6H)-one
15. Cedr-8-ene
16. Murolan-3,9(11)-diene-10-peroxy
17. α -Cadinol
18. Cadalene
19. Coumarin

*According to library search results

Initial use of MonoTrap...

Introduction to the convenient MonoTrap Start UP-KIT

	Description	Qty.
①	MT Holder	5
②	MT Stand	1
③	MT Extract Cup with Vial (20 mL)	5
④	Clean Pin Hole Septum with vial (40mL)	5
⑤	200 μ L glass insert (flat bottom)	40
⑥	MonoTrap DCC18	20
⑦	MonoTrap RCC18	20
⑧	MonoTrap DSC18	20
⑨	MonoTrap RSC18	20



GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome
Shinjuku-ku, Tokyo,
163-1130, Japan
Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences B.V.

De Sleutel 9
5652 AS Eindhoven
The Netherlands
Phone: +31 (0)40 254 95 31
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
Phone: 310-265-4424
Fax: 310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,
Far East International Plaza,
NO.317 Xianxia Road,
Changning District.
Shanghai, China P.C. 200032
Phone: +86 (0)21-6278-2272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn

International Distributors

Visit our Website at www.glsciences.com/distributors